# **Bangladesh University of Engineering and Technology**

Department of Mechanical Engineering



# ME 266 (IPE) Thermal Engineering and Heat Transfer Sessional

Contact Hour: 3.0, Credit Hour: 1.5

# Name of Experiments:

- 1.(a) Determination of Thermal Conductivity of Metals (Copper and Aluminum)
- 1.(b) Study of Thermal Contact Resistance.
- 2. Study of Forced Convection Heat Transfer Over a Flat Plate
- 3. Study of Radiation Heat Transfer

# **Instructions for Students**

# Necessary points of report writing:

- Objectives
- Apparatus (write specification like type, dimensions, range, material, etc.)
- Schematic Diagram of the Experimental Setup (label properly)
- Data and Calculation sheets
- Plotting graphs if any
- Sample Calculation for a particular observation
- Results and Discussions (Result and graph analysis, sources of error, means of improvement, question and answers if any)
- Conclusion

# **Important Instructions:**

- $\checkmark$  Submit the reports in plastic spiral bound files of color specified for the group.
- ✓ You must come to the class before the starting time (for example 2.30 pm). <u>No</u> <u>late attendance.</u>
- $\checkmark$  You must submit all the reports of a group together before the end of a lab session.
- ✓ You must come to the class with prior preparation (study the lab sheet supplied, bring necessary graph papers, pages, calculator, reference books, etc.). <u>You must</u> bring this lab instruction sheet during the sessional class.
- $\checkmark$  Viva on the individual experiments will depend on the concerned teacher.
- ✓ Copying another student's report will be severely dealt with.

# Experiment No. 1 a) Determination of Thermal Conductivity of Metals (Copper and Aluminum) and b) Study of Thermal Contact Resistance.

#### Introduction

When a temperature gradient exists in a body, heat transfer occurs from the high temperature region of the body to its low temperature region. According to Fourier's law of heat conduction, heat transfer through a homogeneous solid body is directly proportional to the area of the section at right angles to the direction of heat flow, and the temperature gradient in that direction. Mathematically, this can be written as,

$$Q \propto A \frac{dT}{dx},$$
 (1)

where,

Q = Heat flow rate by conduction through the material (W)

A = Surface area of the section at right angle to the direction of heat flow  $(m^2)$ 

dx = Differential length of the specimen in the direction of heat flow (m)

dT = Temperature difference in the differential length of the specimen dx (°C or K)

dT/dx = Temperature gradient in the *x*-direction (K/m or °C/m)

The above equation can be written in the following form,

$$Q = -kA\frac{dT}{dx},\tag{2}$$

where, the proportionality constant k is a transport property known as thermal conductivity (W/mK). It is a measure of the ability of a substance to conduct heat and provides the characteristics of the wall material through an indication of the rate at which energy is transferred by diffusion process. The value of thermal conductivity of a specimen depends on the physical structure of matter, atomic and molecular, its composition and temperature. The temperature gradient dT/dx in Eq. (2) is always negative along the direction of heat flow; because heat transfer occurs from the region of high temperature to that of low temperature of the body. Hence, a minus sign is inserted at the right hand side of Eq. (2) in order to make the heat flow rate Q to be positive.

Thermal conductivity is a thermo-physical property of the material, which is, in general, a function of both temperature (*T*) and location (*s*); i.e., k = f(T, s). For isotropic materials, the value of *k* is the same in all directions, i.e., k = f(T). However, for anisotropic materials such as wood, laminated materials, etc., the value of *k* will have a strong directional dependence.

For some materials over certain temperature range, the variation of thermal conductivity with temperature is almost negligible. The development of experimental approximations of boundary value problems is needed to measure the thermal conductivity. Direct measurement of thermal conductivity is based on the widely used 'steady-state method'. It can provide high accuracy and simple data reduction, however, requires a relatively long time to reach steady state.

In thermal engineering, thermal contact resistance represents the resistance to the heat conduction between two solid bodies. When components are bolted or otherwise pressed together, a knowledge of the thermal performance of such joints are also needed. In these composite systems, the temperature drop across the interface between materials may be appreciable

### Objectives

- 1. To plot the measured temperature versus distance curve, and hence, to determine the temperature gradient.
- 2. To determine thermal conductivity of the metal specimen.
- 3. To compare the calculated thermal conductivity of the metal specimen with its standard value.
- 4. To study the significance of thermal contact resistance.

### **Experimental Setup**

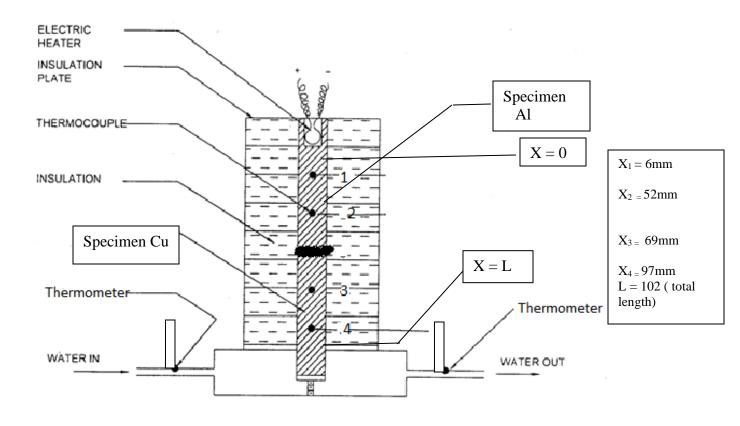


Fig 1. Schematic Diagram of the Experimental Setup.

A schematic diagram of the experimental setup is shown in Fig. 1. The apparatus consists of a vertical stack of specimens clamped between an electrically heated source at the top and a water cooled base, all contained within a Dewar vessel and furnished with a radiation shield and anticonvection baffle. The specimens are fitted with very small thermocouples at known distances apart and connected via a selector switch to a digital temperature readout. The heating current is supplied from a variable voltage power pack and displayed on a digital ammeter. The water cooled base is designed as a calorimeter to measure the heat flow and fitted with very accurate thermometers in the water circuit. Cooling water is supplied from the constant head water tank mounted above the Dewar vessel. A selection of cylindrical specimens of different materials is provided. A Copper specimen is given of 25mm diameter and 38mm long and a Aluminum specimen is given of 25mm diameter and 64 mm long both having cross sectional area of  $4.9 \times 10^{-4} \text{ m}^2$ .

#### **Experimental Procedures**

1. The apparatus has been assembled with one short specimen (Cu) in the lower position and one long specimen (Al) in the upper position.

- 2. The clamping lever being held positioned on the front of the apparatus in the downward position the specimen is placed between the heater and calorimeter block, and clamped in the position by releasing the lever.
- 3. The thermocouple has been inserted in the sequence that they are wired to the support posts, into the holes provided in the sample. The potentiometer-measuring instrument has been connected to the terminals provided on front of the panel.
- 4. The Dewar vessel has been placed on position over the specimens.
- 5. The thermometers has been fitted into the special leak proof connections provided on the top of the calorimeter base (left hand water out, right hand water in), and the water pipes from the water supply to the header tank, the header tank to the inlet on the apparatus, the apparatus outlet connection to drain, via the spring valve provided and the header tank overflow to drain.
- 6. Turn on the water supply and adjust the flow at source to give a small regular overflow from the constant head tank to drain. Adjust the height of the header tank and the clip on the outlet hose to obtain a water flow through the apparatus of 0.5 to 1 cc/sec., whilst maintaining the overflow. During the experiment, if necessary, readjust the clip on the outlet hose to prevent the difference in temperature between the two mercury-in-glass thermometers from exceeding 10°C, whilst maintaining the small overflow to drain.
- 7. The supply voltage has been checked as indicated on the serial number label positioned on the back of the apparatus is correct. The apparatus has already been connected to a single phase AC supply point using the socket provided in the right hand side of the apparatus. The unit is then switched on.
- 8. The heat delivered to the sample is controlled by regulating the current supplied to the heater block using the control knob positioned on the front panel under the ammeter. Turn the knob fully clockwise so that the maximum current is supplied to the heater until temperature T<sub>4</sub>, as indicated by the thermocouple selection knob on the front panel approaches to 80°C and maintained at this temperature until the thermocouples indicate a stable output.

#### **Governing Equations**

The generalized three-dimensional heat conduction equation for a solid body with constant thermal conductivity in Cartesian coordinate can be expressed as:

$$\frac{\partial^2 T}{\partial x^2} + \frac{\partial^2 T}{\partial y^2} + \frac{\partial^2 T}{\partial z^2} + \frac{q'''}{k} = \frac{1}{\alpha} \frac{\partial T}{\partial t},$$
(3)

where,

T = Temperature distribution at any location in the body (°C or K)

x, y, z =Cartesian coordinates (m)

q''' = Internal heat generation rate per unit volume (W/m<sup>3</sup>)

k = Thermal conductivity of the body (W/mK)

 $\alpha$  = Thermal diffusivity [=  $k/\rho C_p$ ] of the body (m<sup>2</sup>/s)

t = Time(s)

 $\rho$  = Density of the body (kg/m<sup>3</sup>)

 $C_p$  = Specific heat at constant pressure (J/kgK)

### **Assumptions**

Following assumptions are made in order to simplify the generalized equation (3) applicable for performing the experiment:

- 1. The heat flow is one-dimensional, i.e., temperature varies along *x*-direction only. This is achieved by putting insulation around the circumferential surface area of the specimen.
- 2. Material is homogeneous and isotropic, and there is no internal heat generation.
- 3. Bounding surfaces of the experimental setup are isothermal in character.
- 4. The temperature gradient in the direction of heat flow is constant, i.e., the temperature profile is linear.
- 5. Heat conduction in steady state condition is achieved before final data is recorded.

### **Derivations of Temperature Distribution**

Under the above assumptions, Eq. (3) can be reduced to the following form,

$$\frac{d^2T}{dx^2} = 0. (4)$$

Integrating Eq. (4) twice with respect to x results,

$$T = c_1 x + c_2, \tag{5}$$

where,  $c_1$  and  $c_2$  are integration constants and can be determined from the appropriate boundary conditions. At steady state condition, the following boundary conditions are satisfied: at x = 0,  $T = T_o$  and at x = L,  $T = T_L$ .

Applying these boundary conditions on Eq. (5), the following expression is obtained,

$$\frac{T - T_o}{T_L - T_o} = \frac{x}{L},\tag{6}$$

where,

T = Temperature of the section at any distance x (°C)

 $T_o$  = Temperature of section at x = 0 (°C)

- $T_L$  = Temperature of section at x = L (°C)
- x = Any distance along the specimen (m)
- L = Distance along the specimen between the sections at x = 0 and x = L (m)

Equation (6) can be rearranged as follows for determination of the theoretical temperature distribution ( $T_t$ ) along the specimen,

$$T_t = \frac{x}{L} \left( T_L - T_o \right) + T_o.$$
<sup>(7)</sup>

#### **Determination of Heat Transfer Rate**

Under steady state condition, the conduction heat transfer along the specimen is equal to the rate of heat carried away by the flowing water. Hence, the heat balance relation becomes,

$$Q_c = Q_w = m_w C_p \Delta T_w, \tag{8}$$

where,

 $Q_c$  = Conduction heat transfer along the specimen (W)

 $Q_w$  = Heat carried away by water (W)

 $m_w$  = Mass flow rate [=  $W_w/t$ ] of water (kg/s)

 $W_w$  = Weight of collected water (kg)

t = time required for water collection (s)

 $C_p$  = Specific heat of water at constant pressure evaluated at  $T_w$  (J/kgK)

 $T_w$  = Mean/bulk temperature [=  $(T_{wo} + T_{wi})/2$ ] of water (°C)

 $\Delta T_w$  = Rise in temperature [=  $T_{wo} - T_{wi}$ ] of flowing water (°C)

 $T_{wi}$  = Water temperature at inlet (°C)

 $T_{wo}$  = Water temperature at outlet (°C)

#### **Determination of Thermal Conductivity**

Using Eqs. (2) and (8), one can obtained the thermal conductivity of a metal as follows:

$$k = -m_{w}C_{p}\Delta T_{w} / A\frac{dT}{dx},$$
(9)

where,

A = Cross-sectional area (=  $\pi D^2/4$ ) of the cylindrical copper specimen (m<sup>2</sup>)

D = Diameter of the cylindrical copper specimen (m)

dT/dx = Temperature gradient obtained from experimentally measured temperature (*T<sub>e</sub>*) versus distance (*x*) curve (W/m)

As per the manufacturing catalogue, the value of thermal conductivity for Aluminum (Al) is 172 W/mk, and for Copper (Cu) is 400 W/mk.

### Discussion

- 1. Briefly explain the nature of experimental temperature distribution plot.
- 2. Is there any deviation in the values of thermal conductivity obtained in the two observations during the experiment? If yes, why?
- 3. Is there any discrepancy between the actual and the experimental values? If yes, why?
- 4. Discuss about thermal contact resistance from the measured temperature vs distance graph.
- 5. What happens to the thermal contact resistance if the spring pressure is increased? Also what will be the scenario if an absolute pressure is given

### Conclusion

Write down the summary of key findings and observations as outlined in the objectives of this experiment.

### **Review Questions:**

- 1. What do you mean by steady state condition and why is it necessary?
- 2. How can you maintain one-dimensional steady state heat conduction?
- 3. How does the value of thermal conductivity of copper vary with the change of its temperature?
- 4. How does the temperature gradient of a solid metal change with its thermal conductivity for a constant heat flux condition?
- 5. Does the conductivity of materials change with the change of water flow rate?

# Data Sheet Determination of Thermal Conductivity of Metals (Cu and Al)

Material of the cylinder is <b>copper and aluminum</b>						
Diameter of the Cu cylinder,	$D_{Cu} =25mm$					
Diameter of the Al cylinder,	$D_{Al} = \dots 25.\dots mm$					

Thermocouple used:					
Туре:	_, Material:	, Range:			

Weight of the empty container, $W_c = \dots$	.kg
Distance between consecutive thermocouple position, $\Delta x = \dots$	.mm
Length of the specimen considered for temperature measurement, $L =102$	mm

# Table 1: Collection of experimental data for different steady state conditions.

		Experimentally measured	Water temp	perature	Weight of container with water, $W_{wc}$ (kg)	Time of collection of water, $t$ (s)
No. of Obs.	Distance, x (m)	temperature (thermocouple reading), $T_e$ (°C)	Inlet T <sub>wi</sub> (°C)	Outlet $T_{wo}$ (°C)		
	1					
1	2					
1	3					
	4					
	1					
2	2					
	3					
	4					

Obs 1 and 2 are for different flow rates of water.

# **Calculation Sheet**

	Obser	rvation
Calculated Parameters	N	lo.
	1	2
$dT/dx$ (°C/m or K/m) [slope of $T_e$ versus x curve]		
Cross-sectional area of the specimen, $A$ (m <sup>2</sup> ) [= $\pi D^2/4$ ]		
Weight of water collected, $W_w$ (kg) [= $W_{wc} - W_c$ ]		
Mass flow rate of water, $m_w$ (kg/s) [= $W_w / t$ ]		
Average water temperature, $T_w$ (°C) [= ( $T_{wo} + T_{wi}$ ) / 2]		
Specific heat of water at $T_w$ , $C_p$ (J/kgK) [use Appendix C]		
Temperature rise of water, $\Delta T_w$ (°C) [= $T_{wo} - T_{wi}$ ]		
Heat carried away by water, $Q_w$ (W) [use Eq.(8)]		
Thermal conductivity of copper, $k_{Cu}$ (W/mK) [use Eq.(9)]		
Thermal conductivity of aluminum, $k_{Al}$ (W/mK) [use Eq.(9)]		
Error (%) [= $ (k_s - k) / k_s  \times 100$ ]		

### Table 2: Calculation of various parameters and thermal conductivity.

Please bring 2 numbers of normal mm graph papers for this experiment.

# **Experiment No. 2** Study of Forced Convection Heat Transfer Over a Flat Plate

#### Introduction

When fluid flows over a solid body or inside a channel, while temperatures of the fluid and the solid surface are different, heat transfer between the fluid and the solid surface takes place as a consequence of the motion of the fluid relative to the surface; this mechanism of heat transfer is called convection. If the fluid motion is artificially induced, say with a pump or a fan that forces the fluid flow over the surface, the heat transfer is said to be by forced convection. If the fluid motion is set up by buoyancy effects resulting from density difference caused by temperature difference in the fluid, the heat transfer is said to be by free (or natural) convection.

There are numerous important engineering applications in which, heat transfer for flow over bodies such as a flat plate, a sphere, a circular tube or a tube bundle are needed.


Depending on the fluid flows over/through the geometry, forced convection can be divided into internal forced convection (flow is confined within a channel, example: flow inside a heated tube) or external forced convection (flow occurs over a surface, example: flow over a flat plate). Fluid flow can also be laminar, transition or turbulent in nature.

When a fluid flows along a surface, irrespective of whether the flow is laminar or turbulent, the particles near the surface are slowed down by virtue of viscous forces. The fluid particles adjacent to the surface stick to it and have zero velocity relative to the boundary (no-slip condition). Other fluid particles attempting to slide over them are retarded because of an interaction between faster and slower moving fluid, a phenomenon that gives rise to shearing forces. The effects of the viscous forces originating at the boundary extend into the body of the fluid, but a short distance from the surface until the velocity of the fluid contained within the region of the substantial velocity change, is called the hydrodynamic boundary layer. The thickness of this boundary layer has been defined as the distance from the surface at which the local velocity reaches 99% of the external velocity (i.e., free steam velocity). The thermal boundary layer also develops along with the hydrodynamic boundary layer due to temperature difference between the surface and the flowing fluid, and the thickness of this boundary layer is defined as the distance from the surface at which the temperature difference between the surface at which the temperature differenc

of the fluid from the surface reaches 99% of the temperature difference between the surface and the free steam value. The relative positions of the hydrodynamic and the thermal boundary layer depend on the value of the Prandtl number. These two boundary layers will coincide with each other, if the value of the Prandtl number is equal to unity. On the other hand, the thermal boundary layer becomes thicker than the hydrodynamic boundary layer when the Prandtl number reaches below unity. Many research have been conducted on forced convection over flat plates and many empirical correlations were proposed depending on the flow and the thermal conditions.

### **Objectives**

- 1. To determine Reynolds number and the critical length of the plate, and to identify the type of flow over the plate.
- 2. To calculate the average convection heat transfer coefficient from thermal balance relation (experimental results) and compare those with the theoretical results (empirical correlations).
- 3. To plot surface temperature distribution along the plate for different heat inputs.
- 4. To plot the variation of average convection heat transfer coefficient (both theoretical and experimental) with the heat input.

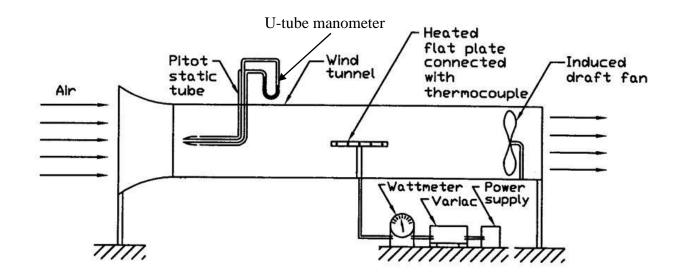

### **Experimental Setup**

Fig 1. Schematic Diagram of the Experimental Setup.

A schematic diagram of the experimental setup is shown in Fig. 1. The main frame of the experimental setup consists of a wind tunnel fitted with an Induced Draft (ID) fan (typeaerofoil, rpm - 2800) to create forced flow through the tunnel, and a horizontal flat plate. The length and the width of the plate are 50.80 and 15.94 cm, respectively and it is placed at the middle of the cross-section of the wind tunnel using a support stand. The flat plate is constructed in such a way that there is a sandwich like heating coil placed inside the plate. As a result, the top and the bottom surfaces of the plate are heated uniformly by the heating coil. Insulation is applied on the surrounding sides of the plate in order to make the heat flow onedimensional.

At the entry of the wind tunnel, a pitot-static tube is placed along the air-flow direction, which is connected to a U-tube manometer filled with water. Heat input is regulated by a variac and is measured by a wattmeter. Thermocouple wires are attached to the heated surface of the plate at five different locations along the plate. Temperature reading from the thermocouples can be controlled using a multipoint selector (rotary type) switch and can be recorded from a digital temperature indicator. Other apparatus of the setup is wall-mounted thermometer to measure surrounding air temperature of the room.

#### **Experimental Procedures**

- 1. Turn on the wind tunnel by switching on the ID fan, and ensure power supply to the heating coil of the flat plate.
- 2. Adjust the variac in order to set the desired power input to the electric heater.
- 3. Observe the temperature readings of the thermocouples and wait for steady state condition. When the steady state is reached, the temperature readings will not change significantly.
- 4. Once the steady state condition is attained, record the temperature reading of the thermocouples from the digital temperature indicator for all five locations of the plate surface.
- 5. Record manometer deflection from U-tube manometer, room temperature from wallmounted thermometer and the supplied power input from wattmeter.
- 6. Repeat steps 2-4 for different power inputs.

#### Assumptions

Following assumptions are considered during the experiment:

1. Steady state is attained before recording the final data.

- 2. Loss of heat by conduction through thermocouple wires and the support stand for the flat plate is neglected.
- 3. Heat loss through the side ends of the plate has been ignored.
- 4. Flow is considered as incompressible and uniform over the flat plate.
- 5. Both hydrodynamic and thermal boundary layers are considered to be developed over the top and the bottom surfaces of the plate thoroughly without being distorted by the support stand.

#### **Governing Equations**

#### Air Velocity in the Wind Tunnel

We know,  $P = \rho g h$ 

Hence, for a certain constant pressure,  $\rho_{air}h_{air} = \rho_{water}h_{water}$ 

$$h_{air} = \frac{\rho_{water} h_{water}}{\rho_{air}},\tag{1}$$

where,

 $h_{air}$  = Equivalent deflection of air in manometer (m)

 $h_{water}$  = Deflection of water in manometer (m)

 $\rho_{air}$  = Density of air at room temperature (kg/m<sup>3</sup>)

 $\rho_{water}$  = Density of water at room temperature (kg/m<sup>3</sup>)

Now, the velocity of air (V) in m/s flowing through the wind tunnel is,

$$V = \sqrt{2gh_{air}}.$$

#### **Governing Parameters for External Forced Convection Heat Transfer**

Table 1: List of governing parameters for external forced convection heat transfer.

Parameters	Definition	Expression	Limiting Value
Reynolds	Inertia Force	$\operatorname{Re}_{L} = \frac{VL}{V}$	$Re_{cr} = 5 \times 10^5$ for flow over a flat
Number	Viscous Force	$KC_L = V$	plate
Prandtl	Momentum Diffusivity	$Pr = \frac{V}{V}$	$Pr \approx 0.7$ for air and
Number	Thermal Diffusivity	$\alpha$	$Pr \approx 4.1 \sim 6.5$ for water (20° ~ 40°C)
Nusselt	Convection	$N\mu = \frac{hL}{hL}$	Nu = 1 for pure conduction
Number	Conduction	$Nu_L = \frac{hL}{k}$	Nu > 1 for convection

Here, v,  $\alpha$ , k are kinematic viscosity (m<sup>2</sup>/s), thermal diffusivity (m<sup>2</sup>/s) and thermal conductivity (W/mK) of fluid, respectively, L is the length of the plate (m) and h is the convection heat transfer coefficient (W/m<sup>2</sup>K),  $Re_{cr}$  is the critical Reynolds number.

Thermo-physical properties of the fluid should be taken at film temperature,

$$T_{f} = \frac{(T_{s} + T_{\infty})}{2}, \qquad T_{s} = \frac{1}{n} \sum_{i=1,2...,n} T_{si}$$
 (3)

where,  $T_s$  is the average surface temperature (°C or K) and  $T_{\infty}$  is the ambient temperature of flowing fluid (°C or K).

#### **Critical Length of the Plate**

It is the length up to which the flow is laminar. The ratio of the critical length  $(x_{cr})$  to the length of the plate (*L*) can be calculated from the following relation:

$$\frac{x_{cr}}{L} = \frac{Re_{cr}}{Re_L} = \frac{5 \times 10^5}{Re_L}.$$
(4)

Using the ratio of  $x_{cr}/L$  and the value of  $Re_L$ , one can identify the type of flow over a flat plate as listed in Table 2.

#### **Empirical Correlations of Average Nusselt Number**

Table 2: List of empirical correlations for forced convection over a flat plate.

Type of	Restrictions	Thermal Boundary Conditions				
flow	Restrictions	Isothermal ( $T_w = \text{constant}$ )	Isoflux ( $q_w = \text{constant}$ )			
Laminar	$Re_{L} < 5 \times 10^{5} \\ 0.6 < Pr < 50 \\ x_{cr}/L > 0.7$	$Nu_L = 0.664 Re_L^{1/2} Pr^{1/3}$	$Nu_L = 0.680 Re_L^{1/2} Pr^{1/3}$			
Turbulent	$5 \times 10^5 \le \text{Re}_L \le 10^7$ $0.6 \le \text{Pr} \le 60$ $x_{cr}/L < 0.3$	$Nu_L = 0.037 Re_L^{4/5} Pr^{1/3}$	$Nu_L = 0.037 Re_L^{4/5} Pr^{1/3}$			
Partly Laminar, Partly Turbulent	$5 \times 10^{5} \le \text{Re}_{\text{L}} \le 10^{7}$ 0.6 \le Pr \le 60 0.3 \le x <sub>cr</sub> /L \le 0.7	$Nu_{L} = (0.037 Re_{L}^{4/5} - 871) Pr^{1/3}$	$Nu_{L} = \frac{0.037 Re_{L}^{4/5} Pr^{1/3}}{1 + 12.33 \times 10^{6} Re_{L}^{-6/5}}$			

Here,  $T_w$  is the wall temperature of the flat plate and  $q_w$  is the wall heat flux applied on the flat plate.

Once the average Nusselt number is calculated from the appropriate correlation listed in Table 2, the <u>theoretical</u> average convection heat transfer coefficient is obtained as follows:

$$h = \frac{Nu_L k}{L}.$$
(5)

#### **Thermal Balance Relation**

The fluid must take away the heat supplied to the plate under steady state condition. Hence, heat balance relation is as follows:

$$Q_s = Q_C + Q_R + Q_K, \tag{6}$$

where,  $Q_s$  – Heat Supplied to the plate (W)

 $Q_C$  – Convection Heat Transfer by the flowing fluid (W)

 $Q_R$  – Radiation Heat Transfer by the flowing fluid (W)

 $Q_K$  – Conduction Heat Transfer through thermocouple wires and support stand for the plate (neglected) (W)

Now, from Newton's law of cooling,

$$Q_C = hA(T_s - T_{\infty}) \tag{7}$$

where,

h = Average convection heat transfer coefficient (W/m<sup>2</sup>K)

A = Total area (=  $2 \times L \times W$ ) of the top and the bottom surfaces of the plate (m<sup>2</sup>)

W = Width of the plate (m)

 $T_s$  = Average surface temperature of the plate (°C or K)

 $T_{\infty}$  = Ambient temperature of the flowing fluid (°C or K)

Similarly, from Stefan-Boltzmann's law,

$$Q_R = \sigma \varepsilon A \left( T_s^4 - T_{\infty}^4 \right), \tag{8}$$

where,

 $T_s$  = Average surface temperature of the plate (K)

 $T_{\infty}$  = Ambient temperature of the flowing fluid (K)

 $\sigma$  = Stefan-Boltzmann constant (= 5.669 × 10<sup>-8</sup> W/m<sup>2</sup>K<sup>4</sup>)

 $\varepsilon$  = Emissivity of the heated plate

Then, using (7) and (8) in Eq.(6), one can get the expression of average convection heat transfer coefficient (<u>experimental</u> result) as follows:

$$h = \frac{Q_s - \varepsilon \sigma A \left(T_s^4 - T_{\infty}^4\right)}{A \left(T_s - T_{\infty}\right)}$$

$$\Rightarrow h = \frac{Q_s - 2LW \varepsilon \sigma \left(T_s^4 - T_{\infty}^4\right)}{2LW \left(T_s - T_{\infty}\right)}.$$
(9)

### Discussions

- 1. Discuss the benefits of calculating average convective heat transfer coefficient over local convective heat transfer coefficient.
- 2. Discuss the nature temperature profile of the flat plate along its length.
- 3. Discuss the effect of variation of power input on average convection heat transfer coefficient.
- 4. Discuss the reasons for difference between the experimental and the theoretical values of the convective heat transfer coefficient.
- 5. Show the development of typical hydrodynamic and thermal boundary layers over the flat plate that you expect from your experimental results.

### Conclusions

Write down the summary of key findings and observations as outlined in the objectives of this experiment.

### **Sample Quiz Questions:**

- 1. What is the function of pitot-static tube?
- 2. What are the advantages of using induced draft (ID) fan over forced draft (FD) fan?
- 3. What will happen when water is flowing over the heated flat plate instead of air? Explain with the concept of hydrodynamic and thermal boundary layers.
- 4. What are the physical significances of Reynolds number, Prandtl number and Nusselt number?
- 5. For forced convection heat transfer, Nu = f(Re, Pr). With this correlation, explain the possible ways to enhance the convection heat transfer coefficient.

# Data Sheet Study of Forced Convection Heat Transfer Over a Flat Plate

Material of the plate is <b>copp</b>	er
Length of copper plate,	$L = \dots $
Width of the copper plate,	$W = \dots $
Thermocouple used:	
Type:, Material:	, Range:
Ambient (room) temperature	e of flowing fluid, $T_{\infty}$ =°C
Density of air at $T_{\infty}$ ,	$\rho_{air}$ = kg/m <sup>3</sup> [use Appendix B]
Density of water at $T_{\infty}$ ,	$\rho_{water} = \dots kg/m^3$ [use Appendix C]
Emissivity of copper plate,	arepsilon = 0.78

Table 1: Collection of experimental data for different observations.

	Power		Thermo	Inlet	Manometer			
No.	Input,	<i>x</i> <sup>1</sup> (m)	<i>x</i> <sub>2</sub> (m)	<i>x</i> <sub>3</sub> (m)	<i>x</i> <sub>4</sub> (m)	<i>x</i> <sub>5</sub> (m)	Air	deflection
of	$Q_{in}$						Temp,	of water,
Obs.	$\mathcal{Q}^{in}$ (W)	No.1	No.2	No.3	No.4	No.5	$T_\infty$	$h_{water}$
	(**)	$T_{s1}$	$T_{s2}$	$T_{s3}$	$T_{s4}$	$T_{s5}$	(°C)	(m)
1								
2								
3								

# **Calculation Sheet**

Table 2: Thermo-physical properties of air at different film temperatures.

Properties	Observation No. 1	Observation No. 2	Observation No. 3
	Film temperature,	Film temperature,	Film temperature,
of air	$T_f = \dots K$	$T_f = \dots K$	$T_f = \dots K$
<i>k</i> (W/mK)			
$\alpha$ (m <sup>2</sup> /s)			
$v (m^2/s)$			

# **Calculated Result**

 Table 3: Calculated governing parameters and convection heat transfer coefficient.

					Average	Convection	n Heat Transfer	
L	<b>_</b> Power Air		Gov	Governing		Coefficient,		
atio.	Input	Velocity	Para	ameter	S		h	
Observation No.	$Q_s$	V				$(W/m^2K)$		
Obs	(W)	(m/s)	$Re_L$	Pr	$\frac{x_{cr}}{L}$	Empirical Co	orrelation	Thermal Balance
					L	Isothermal	Isoflux	Experimental
1								
2								
3								

Please bring two normal mm graph papers for this experiment.

# Appendix A

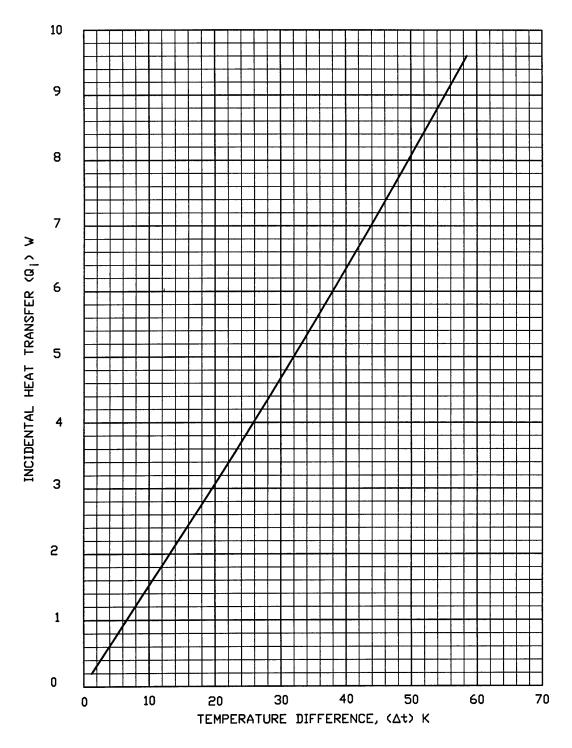


Figure: A. Calibration Curve for Incidental Heat Transfer with Temperature Difference.

# Appendix B

Temperature		TemperatureDensityof therm $\rho$ expansion		Coefficient of thermal expansion	Specific heat $C_p$	Thermal conductivity k	Thermal diffusivity $\alpha \times 10^6$	Absolute viscosity $\mu \times 10^6$	Kinematic viscosity $\nu \times 10^6$
°F	K	°C	kg/m <sup>3</sup>	$\beta \times 10^3$ 1/K	J/kg K	W/m K	m²/s	Ns/m <sup>2</sup>	m²/s
32	273	0	1.252	3.66	1011	0.0237	19.2	17.456	13.9
68	293	20	1.164	3.41	1012	0.0251	22.0	18.240	15.7
104	313	40	1.092	3.19	1014	0.0265	24.8	10.123	17.6
140	333	60	1.025	3.00	1017	0.0279	27.6	10.907	19.4
176	353	80	0.968	2.83	1019	0.0293	30.6	20.790	21.5
212	373	100	0.916	2.68	1022	0.0307	33.6	21.673	23.6
392	473	200	0.723	2.11	1035	0.0370	49.7	25.693	35.5

# Thermodynamic Properties of Dry Air at Atmospheric Pressure

# Appendix C

# Thermodynamic Properties of Water at Saturation Pressure

-			Density	Specific heat	Thermal conductivity	Thermal diffusivity	Absolute viscosity	Kinematic viscosity
T °F K °C		hokg/m <sup>3</sup>	с <sub>р</sub> J/kg К	k W/m K	$lpha  imes 10^6$ m <sup>2</sup> /s	$\mu  imes 10^6$ Ns/m <sup>2</sup>	$v \times 10^6$ m <sup>2</sup> /s	
32	273	0	999.9	4226	0.558	0.131	1794	1.789
41	278	5	1000	4206	0.568	0.135	1535	1.535
50	283	10	999.7	4195	0.577	0.137	1296	1.300
59	288	15	999.1	4187	0.585	0.141	1136	1.146
68	293	20	998.2	4182	0.597	0.143	993	1.006
77	298	25	997.1	4178	0.606	0.146	880.6	0.884
86	303	30	995.7	4176	0.615	0.149	792.4	0.805
95	308	35	994.1	4175	0.624	0.150	719.8	0.725
104	313	40	992.2	4175	0.633	0.151	658.0	0.658
113	318	45	990.2	4176	0.640	0.155	605.1	0.611
122	323	50	988.1	4178	0.646	0.157	555.1	0.556

Example : Here, if  $\mu \times 10^6 = 1.781$ , then  $\mu = 1.781 \times 10^{-6} N s / m^2$ 

Again if  $v \times 10 = 1.785$ , then  $v = 1.785 \times 10^{-1} m^2/s$ 

# **Experiment No. 3 Study of Radiation Heat Transfer**

#### Introduction

Heat transfer from a hot solid body to the surroundings takes place mainly by two processes: convection and radiation. Radiation is the heat transfer by the emission of electromagnetic waves, which carries energy away from the emitting object. For ordinary temperatures (less than red-hot), the radiation is in the infrared region of the electromagnetic spectrum. Heat loss by radiation does not depend on the nature of the surrounding fluid. Radiation may take place through perfect vacuum. However, heat transfer by convection depends on the conditions of the surrounding fluid.

If a hot object (electrically heated element) radiates energy at a temperature  $T_E$  to its cooler surroundings (inside a vessel) at a temperature of  $T_V$ , the net heat transfer by radiation can be calculated with the help of the Stefan-Boltzmann Law as given below:

$$Q_R = \varepsilon A \sigma \left( T_E^4 - T_V^4 \right), \tag{1}$$

where,

 $Q_R$  = Radiation heat transfer from the hot body (W)

 $T_E$  = Absolute temperature of the body (K)

 $T_V$  = Absolute temperature of the surrounding (K)

A = Exposed area of the body (m<sup>2</sup>)

 $\varepsilon$  = Emissivity of the surface of the body

 $\sigma$  = Stefan-Boltzmann constant (= 5.669 × 10<sup>-8</sup> W/m<sup>2</sup>K<sup>4</sup>)

#### Objectives

- 1. To determine temperature difference  $(T_E T_V)$  at zero pressure by plotting  $(T_E T_V)$  versus  $H^{1/4}$  graph.
- 2. To determine the emissivity of the electrically heated element.
- 3. To calculate the radiation heat transfer.
- 4. To analyze the variation of radiation heat transfer with  $(T_E T_V)$ , and hence, to calculate radiation heat transfer coefficient.
- 5. To calculate the heat loss by convection.

#### **Experimental Setup**

A schematic of the experimental setup is shown in Fig. 1. The setup consists of a cylindrical pressure vessel (of diameter 450 mm and height 465 mm) containing a solid cylindrical element (6.35 mm in diameter and 160 mm in length) as a test specimen. The vessel may be charged with air or other gas at wide range of pressures. The element is suspended horizontally from the top cover-plate of the vessel and is finished with a matt black surface. The top cover plate, from which the element is suspended, is fixed with bolts. The solid element is heated internally by means of glass-insulated electric heater and its surface temperature is measured by a thermocouple at a certain point.

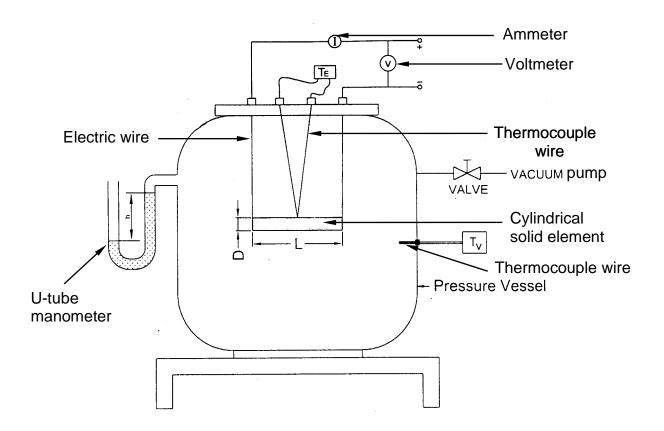


Fig 1. Schematic Diagram of the Experimental Setup.

The solid element is located at sufficiently remote from the walls of the vessel to get substantial free convection. The heat input to the element may range up to about 10W and the maximum working temperature is 200°C. With this small heat input, heating of the pressure vessel walls is considered to be negligible, and the temperature of the surrounding air/gas inside the vessel may be taken as equal to that of the inner wall of the vessel, which is measured by a thermocouple fitted to the vessel wall. The pressure vessel is connected by a

copper tube and an isolating valve to an electrically driven vacuum pump (single stage, rotary type, capacity: 84 liters/min). The pressure in the vessel is determined by the mercury U-tube manometer. Other apparatus of the setup are barometer, voltmeter, ammeter, digital temperature indicator, rheostat, isolating control valve, selector switch, etc.

#### **Experimental Procedures**

- 1. The experiment will start with switching on the vacuum pump to evacuate the vessel until the pressure in the vessel is reduced to the minimum attainable absolute pressure.
- 2. Then the heater will be turned on adjusting the rheostat for a power input of about 5W.
- 3. Next, the isolating valve will be closed, and the vacuum pump will be shut off.
- 4. Now allow sufficient time in order to achieve the steady state condition, which will be indicated from the temperature readings of the thermocouples. When the temperature readings become almost constant with respect to time, the steady state is attained.
- 5. Note down the temperatures of the element and the vessel. Also, record the manometer deflection reading, voltmeter and ammeter readings.
- 6. Repeat the experiment at progressively higher pressures by letting the surrounding air to enter the vessel with the help of slowing opening the isolating control valve. The final reading will be taken at fully open condition of the control valve, when the pressure inside the vessel is equal to the atmospheric pressure.
- 7. In order to maintain constant heat input during each observation, check the voltmeter and the ammeter readings to ensure that they have remained approximately constant. This can be achieved by adjusting the rheostat if necessary.

### **Assumptions**

Following assumptions are considered during the experiment:

- 1. Steady state condition is maintained before recording the final data.
- 2. Readings recorded by voltmeter and ammeter are kept constant throughout the experiment.
- 3. Power loss due to electrical resistance of the leads remains constant at 4 percent.

### **Governing Equations**

### **Heat Supplied to the Element**

Since a power loss of 4% of the product *VI* is considered as a consequence of the electrical resistance of the leads that supply power and support it, the expression of heat input to the solid element becomes,

$$Q_{\rm s} = 0.96VI, \tag{2}$$

where,

 $Q_s$  = Supplied heat input to the element (W)

V =Voltmeter reading (V)

I =Ammeter reading (A)

#### Heat Loss by Conduction

A small part of the heat supplied to the element is lost through conduction along the leads and along the thermocouple wires. This loss is equivalent to the following approximation:

$$Q_{K} = 0.0017 \left( T_{E} - T_{V} \right), \tag{3}$$

where,

 $Q_K$  = Conduction heat transfer along thermocouple wires and along the leads (W)

 $T_E$  = Temperature of the solid cylindrical element (°C or K)

 $T_V$  = Temperature of the vessel wall (°C or K)

#### **Equivalent Surface Area**

Due to the connection of the current carrying leads and the thermocouple leads to the solid cylindrical element, the surface area of the element including the end surfaces may be calculated using the following relation,

$$A = 1.02 \left( \frac{\pi D^2}{2} + \pi DL \right),\tag{4}$$

where,

A = Total exposed surface area of the element  $(m^2)$ 

D = Diameter of the cylindrical solid element (m)

L = Length of the cylindrical solid element (m)

#### Heat Loss by Convection

The surrounding fluid must take away the heat supplied to the element. Hence, heat balance relation is as follows:

$$Q_S = Q_C + Q_R + Q_K, \tag{5}$$

where,  $Q_C$  is convection heat loss by the surrounding fluid (W). Now, using (1), (2) and (3) in Eq.(5), one can get the expression of convection heat loss as follows:

$$Q_{c} = Q_{s} - Q_{R} - Q_{K} = 0.96VI - \varepsilon A\sigma \left(T_{E}^{4} - T_{V}^{4}\right) - 0.0017 \left(T_{E} - T_{V}\right).$$
(6)

#### **Emissivity of the Element**

If the pressure vessel can be maintained in perfect vacuum condition, there will be no heat loss by convection, and then Eq.(6) is reduced to as follows:

$$0 = 0.96VI - \varepsilon A\sigma \left(T_{E}^{4} - T_{V}^{4}\right) - 0.0017 \left(T_{E} - T_{V}\right)$$
  
$$\Rightarrow \varepsilon = \frac{0.96VI - 0.0017 \left(T_{E} - T_{V}\right)}{A\sigma \left(T_{E}^{4} - T_{V}^{4}\right)}.$$
(7)

In order to determine the emissivity of the element ( $\varepsilon$ ) from Eq.(7), it is necessary to measure temperatures of the element and the vessel at zero pressure (perfect vacuum condition). However, due to limitation of experimental setup, it is not possible to attain perfect vacuum condition. Hence, an alternate procedure is followed for calculating the emissivity as mentioned in the following steps:

- 1. Plot temperature difference  $(T_E T_V)$  versus  $H^{1/4}$  graph (where, H is the absolute pressure in mmHg) and determine  $(T_E T_V)$  at zero pressure from extrapolation.
- 2. Assume  $T_V$  at the lowest pressure reading as the value of  $T_V$  at zero pressure. Hence, calculate  $T_E$  at zero pressure from the extrapolation data of  $(T_E T_V)$  at zero pressure as obtained in the previous step.
- 3. Finally, insert the values of  $T_E$ ,  $T_V$ , V, I,  $\sigma$ , and A from Eq.(4) in Eq.(7) to calculate the emissivity of the element.

#### **Radiation Heat Transfer Coefficient**

Since  $Q_R$  versus  $(T_E - T_V)$  graph is almost linear, one can consider the following relationship analogous to Newton's law of cooling for convection heat transfer:

$$Q_R = h_R A \left( T_E - T_V \right), \tag{8}$$

where,  $h_R$  is the radiation heat transfer coefficient (W/m<sup>2</sup>K) and can be obtained from the slope of the  $Q_R$  versus ( $T_E - T_V$ ) curve as follows:

$$h_R = \frac{slope}{A}.$$
(9)

#### Discussions

- 1. Briefly explain the nature of  $(T_E T_V)$  versus  $H^{1/4}$  graph.
- 2. Discuss the effects of temperature difference on radiation heat transfer from  $Q_R$  versus  $(T_E T_V)$  graph obtained from the experiment.
- 3. Discuss on the value of radiation heat transfer coefficient  $(h_R)$  obtained from the experiment.

4. Discuss the reasons for difference between the expected value and the calculated value of emissivity obtained from the experiment.

### Conclusions

Write down the summary of key findings and observations as outlined in the objectives of this experiment.

### **Sample Quiz Questions:**

- 1. What is emissivity and how does its value vary from black body to white body?
- 2. What are the limitations of calculating emissivity of the element observed during the experiment?
- 3. Mention the factors on which the rate of emission of radiation depends.

# Data Sheet Study of Radiation Heat Transfer

Diameter of the solid element,	D =m	
Length of the solid element,	L =m	
Thermocouple used:		
Type:, Material:	, Range:	

Barometric Pressure, P<sub>atm</sub> = .....mm Hg

	Supply Supply U-tube Manomete		U-tube Manometer	Element	Vessel	
No. of	Voltage,	Current,	reading,	temperature,	temperature,	
Observation	V	Ι	h	$T_E$	$T_V$	
	(V)	(A)	(mmHg)	(°C)	(°C)	
1						
2						
3						
4						
5						
6						
7						
8						

# Table 1: Collection of experimental data for different observations.

# **Calculation Sheet**

Coloulated Devenuetors	Number of Observations								
Calculated Parameters	1	2	3	4	5	6	7	8	
$T_E(\mathbf{K})$									
$T_V(\mathbf{K})$									
$T_E - T_V(\mathbf{K})$									
$H = P_{atm} - h \text{ (mm Hg)}$									
$H^{1/4} (\mathrm{mm \ Hg})^{1/4}$									
$A (m^2)$ [use Eq.(4)]									
$\varepsilon$ [use Eq.(7)]									
$Q_{S}(W)$ [use Eq.(2)]									
$Q_K$ (W) [use Eq.(3)]									
$Q_R$ (W) [use Eq.(1)]									
$Q_C$ (W) [use Eq.(6)]									
$h_R$ (W/m <sup>2</sup> K) [use Eq.(9)]									

 Table 2: Calculated various parameters and emissivity of the element.

Please bring two mm graph papers for this experiment.